


RockImager: Putting a plate in and removing it

RockImager software controls the RockImagers. And it is installed on two computers. Each install controls either the 20 °C RockImager or the 8 °C RockImager.

Users will have little need to use the RockImager software except when entering and retrieving an experimental plate.


The RockImager software should never be closed.

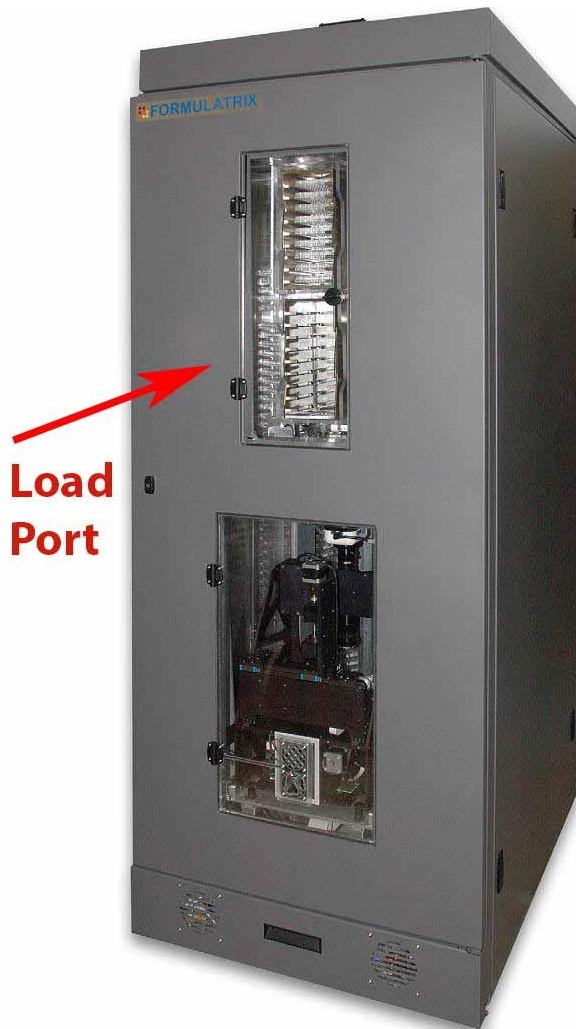

Putting an experiment plate in the Rock imager

Open the door of the Load Port and place your plate in the hotel with the barcode facing inwards. Close the door and go to the RockImager software on the computer next to the RockImager you are using.

From the Load Port menu select Store Contents. The RockImager will scan the Load Port hotel for plates and display their bar coded information in the Load Port tab on the Plates tab. When this information is displayed for your plates you can walk away, the plates will be put in storage and imaged according to the schedule you set.

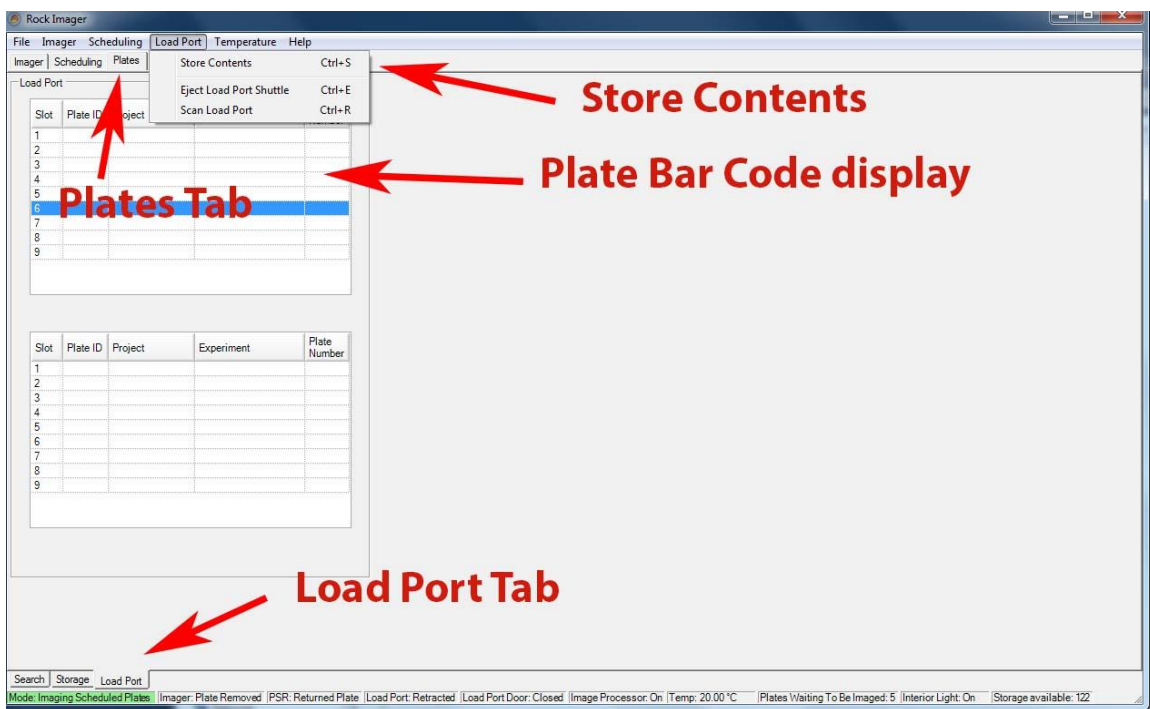

Searching for a plate you have stored in a RockImager

The RockImager software allows you to find your plate in the RockImager. From the Plates Tab select the Search Tab. The top line (in blue) is where you put the search target each your Plate ID. When your plates are found they will be displayed in the search field.

